

*Simultaneous isolation and concentration with SmartFlow™ TFF*


Each parameter of the TFF process: product, membrane type, shear, pore size, temperature, concentration factor, pH, anti-foam, etc. may impact the fermentation broth components passage through the membrane. This is why a systematic experimental plan must be developed and executed to optimize a concentration and diafiltration process.

- 1) Select the *SmartFlow™* filter modules to evaluate for the isolation and concentration steps. The selection requires specifying a combination of membrane type, channel height, and membrane area for a given module that will be tested.
  - a) NCSRT has filtered thousands of solutions and therefore can provide several membrane chemistries and pore sizes that will likely work in the majority of cases. In general the pore size should be 5 to 10 times the size of the molecule to be passed through the membrane and one half to one third the size of a molecule to be retained.
  - b) The combination of the channel height and the fluid velocity through the flow channel created by the recirculation pump produce a shear at the membrane surface. It is this shear that governs the separation performance and efficiency in ultrafiltration steps. Velocity governs the performance in microfiltration steps. Care must be taken in selecting and maintaining the shear and the velocity at the membrane surface.
  - c) The membrane area also affects the pump size required to achieve the necessary shear rates for a given separation.
- 2) Select the first membrane to test for the isolation or cell harvest step. Use Table 1 to determine the recommended starting membrane pore size to pass the desired product. For more specific recommendations, see the specialized Optimization Procedure for your class of target molecule.

Table 1. Suggested membranes by target molecular weight

| Target Molecule MW | Membrane polymer | Pore size |
|--------------------|------------------|-----------|
| < 5 kD             | RC               | 30 kD     |
| 5-30kD             | RC               | 100 kD    |
| 30 - 100 kD        | MPS              | 0.1 um    |
| > 100 kD           | MPS              | 0.2 um    |
| Mab                | MPS              | 0.45 um   |
| < 5 kD             | PES              | 50 kD     |
| 5-10 kD            | PES              | 100 kD    |

- 3) Select the channel height for the module.
  - a) For the isolation of a secreted protein from microbial cells, a channel height between 0.75 and 1.5 mm is recommended.
  - b) In most cases, a channel height of 0.75 mm is recommended because it will require the lowest recirculation (and thus the smallest pump) and produce the highest flux rate.
  - c) Cases to use a higher channel height:
    - i) If cell aggregation is occurring, the lower height channels may clog. If the channel is clogged by aggregates or process particles, the inlet pressure will increase dramatically and the permeate rate will decrease over a short period of time. This will occur usually in the first five minutes.

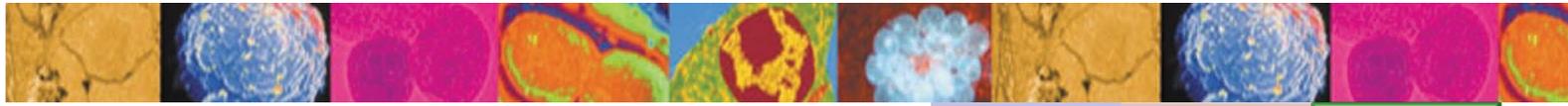

Works™  
Protocol

Works™  
Case Study

Works™  
Optimization  
Procedure

### Simultaneous isolation and concentration with SmartFlow™ TFF

- ii) In cases where high solids are desired, a channel height of 0.875 mm or above will be necessary.
  - 4) Select the membrane area.
    - a) The membrane area depends upon the batch size to be processed. For filtration process development trials, usually the smallest size membrane and thus the smallest batch size is desired.
    - b) For cell harvests an important parameter is the membrane capacity or LM ratio. The membrane capacity or LM ratio is defined as the volume of starting material divided by the membrane area.
    - c) The range of the LM ratio for the isolation of proteins by continuous diafiltration is between 25 to 120 LM.
      - i) The typical starting ratio is 60 LM. See below for additional information on optimizing the LM ratio.
      - ii) If a fermentation broth is being concentrated, the presence of antifoam is an important parameter that can impact the starting volume to membrane area ratio. Antifoam agents may significantly reduce the permeate flux observed with ultrafiltration membranes. In the cases of fermentation broths containing antifoam, use a starting ratio of 30 LM.
    - d) The minimum batch size is the system hold up volume times the concentration factor. For a continuous diafiltration, the minimum batch size is simply the hold up volume.
    - e) The membrane area needed is the batch size divided by the LM ratio.
  - 5) Determine the velocity and shear rate.
    - a) Isolation steps should be optimized on the fluid velocity and concentration steps should be optimized on shear.
    - b) The typical velocity for the isolation of products using microfiltration ranges from 50 cm/sec to 150 cm/sec.
    - c) The benefit of increasing the velocity in an isolation step is an increased passage rate.
    - d) The disadvantages of increasing the channel velocity in a microfiltration are:
      - i) Higher pump costs due to higher recirculation flow rates.
      - ii) Higher pressure drops and TMPs which may decrease the passage of the desired protein.
      - iii) Increased channel velocity may be detrimental to shear sensitive cells such as insect or other mammalian cell lines.
    - e) An increase in the flow rate should be balanced by an increase in the protein passage for the process to retain the same overall efficiency. The energy costs of running the pump at a higher shear rate must be offset by savings on membranes to make increasing the flow rate efficient.
  - 6) Calculate the flow rate needed to operate the selected module at the selected shear rate or velocity using the *WORKS™* Scale-UP LPM GPM spreadsheet. Ensure that a pump is available that can produce this flow rate the needed pressure. If a suitable size pump is not available, consider either running a smaller trial or calling NCSRT to determine if a suitable size pump is available.
  - 7) Use Table 2 to determine the module(s) part numbers for ordering.
  - 8) Select the first membrane to test for the concentration or UF step. Use Table 3 to determine the recommended starting membrane pore size to retain the desired product. For more specific recommendations, see the specialized optimization procedure for your molecule.

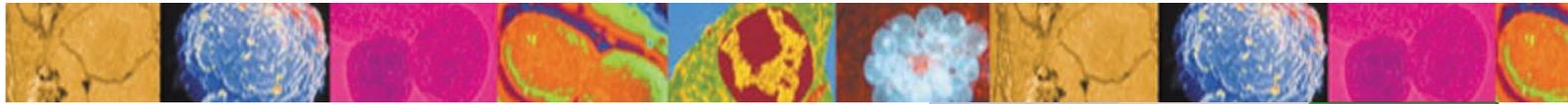



Table 2: SmartFlow™ filter module part numbers

| Module Size                           | Channel Height | Membrane polymer and pore size |
|---------------------------------------|----------------|--------------------------------|
| 74 100 ft <sup>2</sup> Optisep® 11000 | D 0.5 mm       | 5B-0030 RC 30 kD               |
| 72 50 ft <sup>2</sup> Optisep 11000   | E 0.75 mm      | 5B-0100 RC 100 kD              |
| 71 10 ft <sup>2</sup> Optisep 11000   | G 0.875 mm     | 1N-9010 MPS 0.1 μm             |
| 41 10ft <sup>2</sup> Optisep 7000     | H 1 mm         | 1N-9020 MPS 0.2 μm             |
| 40 5 ft <sup>2</sup> Optisep 7000     | J 1.5 mm       | 1N-9045 MPS 0.45 μm            |
| 52 2 ft <sup>2</sup> Optisep 3000     |                | 24-0050 PES 50 kD              |
| 51 1 ft <sup>2</sup> Optisep 3000     |                | 1B-0100 PS 100 kD              |

Table 3. Suggested membranes by target molecular weight

| Target molecule | Membrane polymer | Pore size |
|-----------------|------------------|-----------|
| > 8 kD          | RC               | 5 kD      |
| > 18 kD         | RC               | 10 kD     |
| >50 kD          | RC               | 30 kD     |
| > 180 kD        | RC               | 100 kD    |
| 3- 8 kD         | TFM              | 1 kD      |
| >7.5 kD         | PES              | 5 kD      |
| 18- 100 kD      | PES              | 10 kD     |
| >100 kD         | PES              | 50 kD     |

- 9) Select the channel height for the module.
  - a) For the concentration of a prefiltered protein, a channel height between 0.75 and 1.5 mm is common.
  - b) In most cases, a channel height of 0.75 mm is recommended because it will require the lowest recirculation (and thus the smallest pump) and produce the highest flux rate.
  - c) In cases where high solids are desired, a channel height of 0.875 mm or above will be necessary.
- 10) Select the membrane area.
  - a) The key to sizing the concentration step is to ensure that the permeate rate is equal in the two steps. Usually this requires sizing providing twice the membrane area for the concentration step than was used in the isolation step.
- 11) Determine the shear rate.
  - a) The shear rate for the isolation of secreted proteins by continuous diafiltration ranges from 7,500 sec<sup>-1</sup> to 20,000 sec<sup>-1</sup>.
  - b) The typical starting shear rate for a process development run is 10,000 sec<sup>-1</sup>.
  - c) The benefit of increasing the shear rate is an increased permeate rate.
  - d) The disadvantage of increasing the shear rate is the higher pump costs due to higher recirculation flow rates.
  - e) An increase in the shear rate should be balanced by an increase in the flux rate for the process to retain the same overall efficiency. The energy costs of running the pump at a higher shear rate must be offset by savings on membranes to make increasing the shear rate efficient.
- 12) Calculate the flow rate needed operate the selected module at the selected shear rate using the *WORKS*™ Scale-UP LPM GPM spreadsheet. Ensure that a pump is available that can produce this flow rate at the

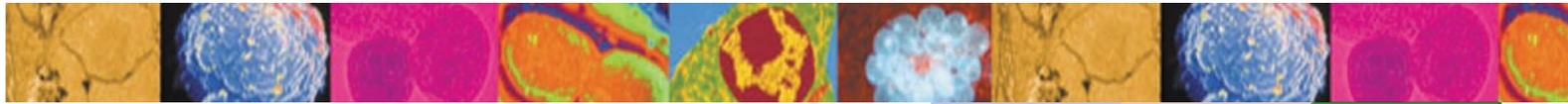

Simultaneous isolation and concentration with SmartFlow™ TFF

Table 4: SmartFlow™ filter module part numbers - concentration step

| Module Size                           | Channel Height | Membrane polymer and pore size |
|---------------------------------------|----------------|--------------------------------|
| 74 100 ft <sup>2</sup> Optisep® 11000 | D              | 0.5 mm 5B-0030 RC 30 kD        |
| 72 50 ft <sup>2</sup> Optisep 11000   | E              | 0.75 mm 5B-0100 RC 100 kD      |
| 71 10 ft <sup>2</sup> Optisep 11000   | G              | 0.875 mm FD-0001 TFM 1 kD      |
| 41 10ft <sup>2</sup> Optisep 7000     | H              | 1 mm 5B-0005 RC 5 kD           |
| 40 5 ft <sup>2</sup> Optisep 7000     | J              | 1.5 mm 5B-0010 RC 10 kD        |
| 52 2 ft <sup>2</sup> Optisep 3000     |                | 2B-0050 PES 50 kD              |
| 51 1 ft <sup>2</sup> Optisep 3000     |                | 2B-0005 PES 5 kD               |

needed pressure. If a suitable size pump is not available, consider either running a smaller trial or calling NCSRT to determine if a suitable size pump is available.

- 13) Use Table 4 to determine the module(s) part numbers for ordering.


Filter Operation:

- 1) After loading the filter modules and making all the connections, the first step is to perform a water and/or buffer rinse of the system directing the permeate to the waste.
- 2) After the rinse, direct the permeate line back to the retentate tank so no concentration occurs prior to establishing the desired shear rate and performing the transmembrane pressure (TMP) optimization procedure.

**IMPORTANT: Do not permit the permeate line to come in contact with the retentate fluid. This can contaminate the permeate pool in later samples.**

- 3) Slowly increase the flow rate recirculation pump to the calculated rate from step 7 above.
- 4) Start with the backpressure at zero. The inlet pressure should be at least 5 psig (0.3 bar) and remain below 12 psig (0.8 bar). If the inlet pressure is above 12 psig (0.8 bar), the recirculation flow rate should be reduced such that the inlet pressure remains below 12 psig (0.8 bar). If the inlet pressure is less than 5 psig, then slight backpressure can be added until the inlet pressure increases to 5 psig. Table 5 provides typical TMP values for the different membranes used in SmartFlow™ TFF filter modules.
- 5) Begin Concentrating.
  - a) Remove the permeate lines from the retentate tank and place them in the permeate vessel.
  - b) Wait until about 5% of the starting retentate volume has passed through the membrane to the permeate to take the initial samples of the retentate and permeate.
  - c) Take permeate and retentate samples when each additional concentration factor is reached.
  - d) When each sample is taken, record the permeate flow rate using a graduated cylinder, scale, or flow meter.
- 6) Once the desired concentration factor is reached in the isolation process, record the volume remaining in the retentate of the isolation loop.
  - a) The remaining volume in the retentate can be calculated by subtracting the permeate volume and retentate volume samples from the starting volumes.

$$\text{Retentate Volume} = \text{Starting Volume} - \text{Permeate Volume} - \text{Retentate Sample Volume.}$$

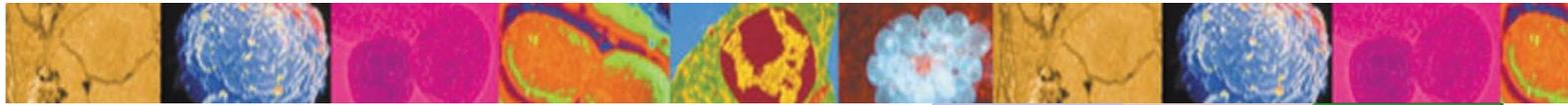


### *Simultaneous isolation and concentration with SmartFlow™ TFF*

- 7) At this point, the concentration skid recirculation pump should be turned on to provide permeate which will act as the buffer for the diafiltration of the isolation step.
  - a) Slowly increase the pump speed until the flow rate calculated in step 12 above is reached.
  - b) Place the permeate hose in the retentate vessel from the isolation step. Here, the permeate will provide the diafiltration buffer for the first step.
  - c) Increase the TMP until it is in the midpoint of the range given in Table 5.
  - d) Take a permeate flow rate from the concentration step permeate.
  - e) If the concentration step permeate is higher than the isolation step permeate, then decrease the pressure until they are equal.
  - f) If the isolation step permeate is higher than the concentration step permeate, then increase the pressure and measure the permeate flow rate.
    - i) Keep increasing the pressure as long as the permeate flow rate is increasing or until it is equal to the isolation step permeate flow rate.
    - ii) If an increase in the pressure does not increase the permeate flow rate (an indication that the gel layer has compressed on the membrane), increase the recirculation rate until the permeate flow rates from the two steps are equal.
- 8) Diafiltration – the following describes the procedure for diafiltering the product 5x:
  - a) Start to monitor the retentate volume in the isolation step tank.
  - b) The level in the tank will remain constant if the two permeate flow rates are equal.
  - c) Continue to add permeate back to the tank until 5 times the retentate volume has been added to the system.
  - d) Take samples from the permeate hose and both the isolation step and concentration step retentate tanks when each diafiltration factor is reached (i.e. take a sample when the permeate volume is equal to a multiple of the retentate volume such as 1X, 2X, etc.).
  - e) For other diafiltration factors, continue the process to the amount of diafiltration buffer equals number of desired diafiltration factor times the system volume recorded in step 6.
    - i) The theoretical recovery from a 3X diafiltration for a molecule with a 100% passage is 95%.
    - ii) Increasing the diafiltration factor will increase the yield especially when the target molecule has low passage. However, the cost of increasing the diafiltration volume is that the process time will be greater and a larger supply of buffer will be needed.
    - iii) Decreasing the diafiltration factor will decrease the yield. However, for molecules with high passage and low value, the small decrease in the yield may be worth the faster processing time and the saving on buffer.

Data analysis:

Sample Analysis:

- 1) Check the permeate samples from the isolation step for cells.
- 2) Use a gel or specific protein assay to check the permeate samples from the concentration step for product passage.
- 3) Calculate the membrane flux rate or LMH (L/m<sup>2</sup>/h) by dividing the measured permeate flow rate at each sample by the membrane area.
- 4) Calculate the instantaneous product passage in the isolation step by dividing the permeate product content by the retentate product content and multiplying by 100.
- 5) Record the data on the Membrane Test worksheet.

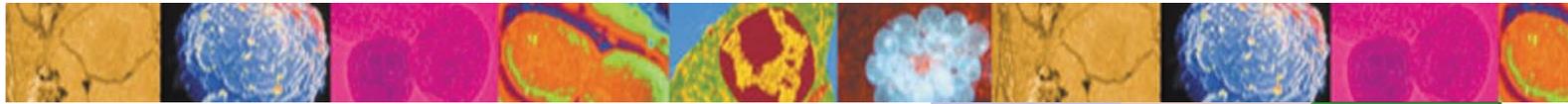

*Simultaneous isolation and concentration with SmartFlow™ TFF*

Process Optimization:


The procedure should be repeated under different process conditions to ensure that the optimized conditions are reached.

- 1) The important variables to optimize are the yield, membrane passage, and membrane flux rate.
- 2) An important parameter that affects the yield, passage, and flux rate for cell harvest is the membrane capacity or LM ratio for the isolation step.
- 3) Increasing the LM ratio decreases membrane performance, which increases processing time and decreases membrane costs. If membrane performance suffers greatly, then saving a little bit on membrane will not offset the costs in higher processing time.
- 4) Decreasing the LM ratio increases the membrane performance and increases membrane costs. Increasing membrane performance may decrease the processing time at a small incremental membrane cost, therefore decreasing total cost.

Table 5 Typical transmembrane pressure values for *SmartFlow* modules

| Membrane Pore Size               | Transmembrane Pressure Starting Value PSIG (Bar) | Transmembrane Pressure Ranges PSIG (Bar) | Cell Harvest Inlet PSIG (Bar) Starting Value | Cell Harvest Outlet PSIG (Bar) |
|----------------------------------|--|--|--|--------------------------------|
| <b>Ultrafiltration Membranes</b> |  |  |  |                                |
| 1 kDa                            | 75 (5)   | 90 to 150 (6 to 10)                      |  |                                |
| 5 kDa                            | 45 (3)   | 60 to 90                                 |  |                                |
| 10 kDa                           | 30 (2)   | 45 to 90 (3 to 6)                        |  |                                |
| 30 kDa                           | 15 (1)   | 30 to 75 (2 to 5)                        |  |                                |
| 100 kDa                          | 15 (1)   | 20 to 60 (1.37 to 4)                     | 20 (1.37)                                    | 12 (0.83)                      |
| 300 kDa                          | 10 (0.69)  | 15 to 45 (1 to 3)                        | 20 (1.37)                                    | 10 (0.69)                      |
| 500 kDa                          | 7.5 (0.5)  | 10 to 30 (0.7 to 2)                      | 7.5 (0.5)                                    | 0                              |
| <b>Microfiltration Membranes</b> |  |  |  |                                |
| 0.1µ                             | 2 (0.13)   | 4 to 15 (0.27 to 1.0)                    | 4 (0.275)                                    | 0                              |
| 0.2µ                             | 2 (0.13)   | 4 to 15 (0.27 to 1.0)                    | 4 (0.275)                                    | 0                              |
| 0.45µ                            | 2 (0.13)   | 4 to 10 (0.27 to 0.69)                   | 4 (0.275)                                    | 0                              |
| 0.8µ                             | 1 (0.07)   | 1 to 6 (0.07 to 0.41)                    | 2 (0.13)                                     | 0                              |
| 1.0µ                             | 1 (0.07)   | 1 to 6 (0.07 to 0.41)                    | 2 (0.13)                                     | 0                              |
| 2.0µ                             | 1 (0.07)   | 1 to 6 (0.07 to 0.41)                    | 2 (0.13)                                     | 0                              |
| 3.0µ                             | 1 (0.07)   | 1 to 6 (0.07 to 0.41)                    | 2 (0.13)                                     | 0                              |

- 5) To find the optimal LM ratio:
  - a) If the current trial was too fast with very high yield, increase the LM ratio by starting with a larger volume of starting material.
  - b) If the current trial was too slow or had a low yield, decrease the LM ratio by starting with a smaller volume of starting material.
- 6) The module used is an important optimization parameter. By changing the membrane chemistry or membrane type, optimized flux rates and passage may be found.

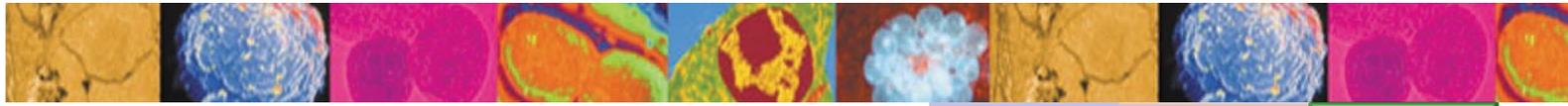

Works™  
Protocol

Works™  
Case Study

Works™  
Optimization  
Procedure

### *Simultaneous isolation and concentration with SmartFlow™ TFF*

- 7) Using the same membrane, the shear rate can be optimized by increasing and decreasing the shear rate and measuring the effects on the membrane flux rate and passage. If an increase in the shear rate results in a relatively large increase in the flux rate, then the savings in membrane cost will offset the increased energy consumption. Likewise, the velocity can be optimized in microfiltration steps. Changes in velocity may affect the passage of the target molecule.
- 8) The concentration factor before starting diafiltration should also be optimized.
- 9) The goal is to begin the diafiltration with the instantaneous passage of at least 50% to increase diafiltration efficiency.
- 10) After analysis of the data, select the best performing modules. The best performing modules will retain the cells, permit the desired media component to pass into the permeate with a high permeate flux for the isolation step. Additionally, the best performing concentration step module will concentrate the desired protein without losses into the permeate.

### Conclusion:

This *SmartFlow™* filter *WORKS™* *Simultaneous isolation and concentration* Optimization Procedure provides guideline for optimizing the application of NCSRT's *SmartFlow* filters. To see how others have applied the technology to their separations operations, please refer to the *Simultaneous isolation and concentration WORKS* Case Study. NCSRT has developed a simple step-by-step protocol that has been proven to deliver >90% product yield in the applications to which it has been applied. Along with the *Simultaneous isolation and concentration WORKS* Protocol, NCSRT has developed a Scale Up Component listing providing the part numbers and ordering information for the *SmartFlow* filter modules to execute the protocol at all process stages and volumes. To receive the complete application package, please request the *Simultaneous isolation and concentration WORK*book.

NCSRT's *SmartFlow* filter technology...It *WORKS*.



NCSRT, Inc.  
1000 Goodworth Dr.  
Apex, NC 27539  
USA

Phone: +1 919 387 8460  
Fax: +1 919 387 8540  
email: [info@ncsrt.com](mailto:info@ncsrt.com)  
Web: <http://www.ncsrt.com>